

FALSE POSITIVES

with BLUESTAR IDENTI-HEM and BLUESTAR OBTI

Tests with beech marten, pine marten, and European polecat blood were carried out.

1) With BLUESTAR IDENTI-HEM

Only EUROPEAN MARTEN blood yielded a slight FALSE POSITIVE.

The test line is present but faint compared to the control line.

The beech marten and pine marten blood did not produce any reaction.



2) With BLUESTAR OBTI

Only EUROPEAN POLECAT blood yielded a FALSE POSITIVE. The test line is similar to the control line.

The beech marten and pine marten blood did not produce any reaction.

