

## FALSE POSITIVE

### with BLUESTAR® IDENTI-HEM® and BLUESTAR® OBTI

---

Tests were carried out using mustelid blood: beech marten, pine marten and European polecat.

#### With BLUESTAR® IDENTI-HEM®

Only European polecat blood caused a light FALSE POSITIVE: The test line (T) is present but very weakly.

Weasel and marten produced no reaction.



#### With BLUESTAR® OBTI

Only European polecat blood produced a FALSE POSITIVE: The test line (T) is similar to the control line (C).

Weasel and marten produced no reaction.

