

Using BLUESPRAY[®] with BLUESTAR[®] Magnum

GETTING READY FOR USE

Unscrew the empty bottle provided in the BLUESPRAY[®] kit.



In its place, screw in the bottle of BLUESTAR[®] Magnum already activated with its oxidant tablets.



SPRAYING

To spray BLUESTAR[®] Magnum, press the push button.



Best results are obtained when releasing pressure **INTERMITTENTLY AND WITHOUT SHAKING** at a temperature between 15 and 40°C (between 60 and 120°F).

If the pressure reservoir gets too cold while holding it in your hand, you reservoir insert it into the supplied foam sleeve which will act as a thermal insulator.

To maintain sufficient pressure during prolonged use, use two pressure reservoirs alternately to replace a cold unit by a warm one.

REPLACING THE RESERVOIR



To replace the pressure reservoir, unlock the used pressure reservoir by turning it counter-clockwise.



Lock in the new pressure reservoir by turning it clockwise on the BLUESPRAY[®] head, while fitting the pushbutton over the valve.

PRECAUTIONS

After use and when travelling, empty or close the BLUESTAR[®] Magnum bottle using the supplied cap.

During handling, do not turn the BLUESPRAY[®] upside down so as to avoid leakage.